



# **Automated Nucleic Acid Isolation from Formalin-Fixed Paraffin-Embedded Tissue and Detection of Cancer Biomarkers**

**Dr. Guido Hennig  
Siemens Healthcare Diagnostics  
Molecular Research Germany**

**Copyright © Siemens Healthcare Diagnostics 2009. All rights reserved.**

# Siemens Healthcare Diagnostics

Molecular Research Germany

**SIEMENS**

## Location

- BioCampus Cologne

## Team of 35 people

- Biologists, chemists, physicians, mathematicians, computer scientists, technicians

## Research Focus


- Molecular (RNA+DNA) diagnostics in cancer
- Nucleic Acid Isolation Technology



## In-Vitro Diagnostic to Guide Breast Cancer Therapy

**Today:** Therapy guidelines based on histopathological parameters leaving still numerous therapy options  
Some molecular testing recently included into guidelines


**Future:** Develop and validate more/better molecular tests to pick optimal therapy guided by molecular profiling of tumor

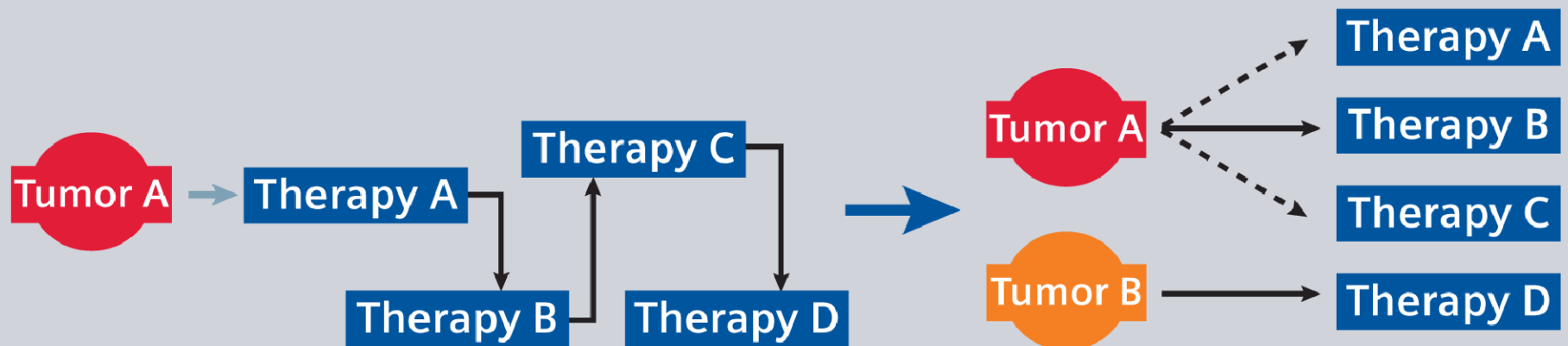

## Why FFPE Tissue for Cancer Diagnostics?

Two principle sources of tumor material exist:


- fresh frozen tissue
- formalin-fixed paraffin-eMBEDDED (FFPE) tissue

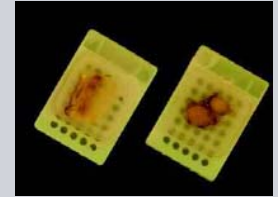


- Access to pathological FFPE archives with linked clinical records allows high-throughput retrospective finding + validation studies
- FFPE tissue is the routine material after tumor surgery and key for successful implementation and commercialization of a diagnostic kit (requires no change of clinical practice)

# Prognostic / Predictive Markers in Breast Cancer

## Research Strategy of Siemens Diagnostics

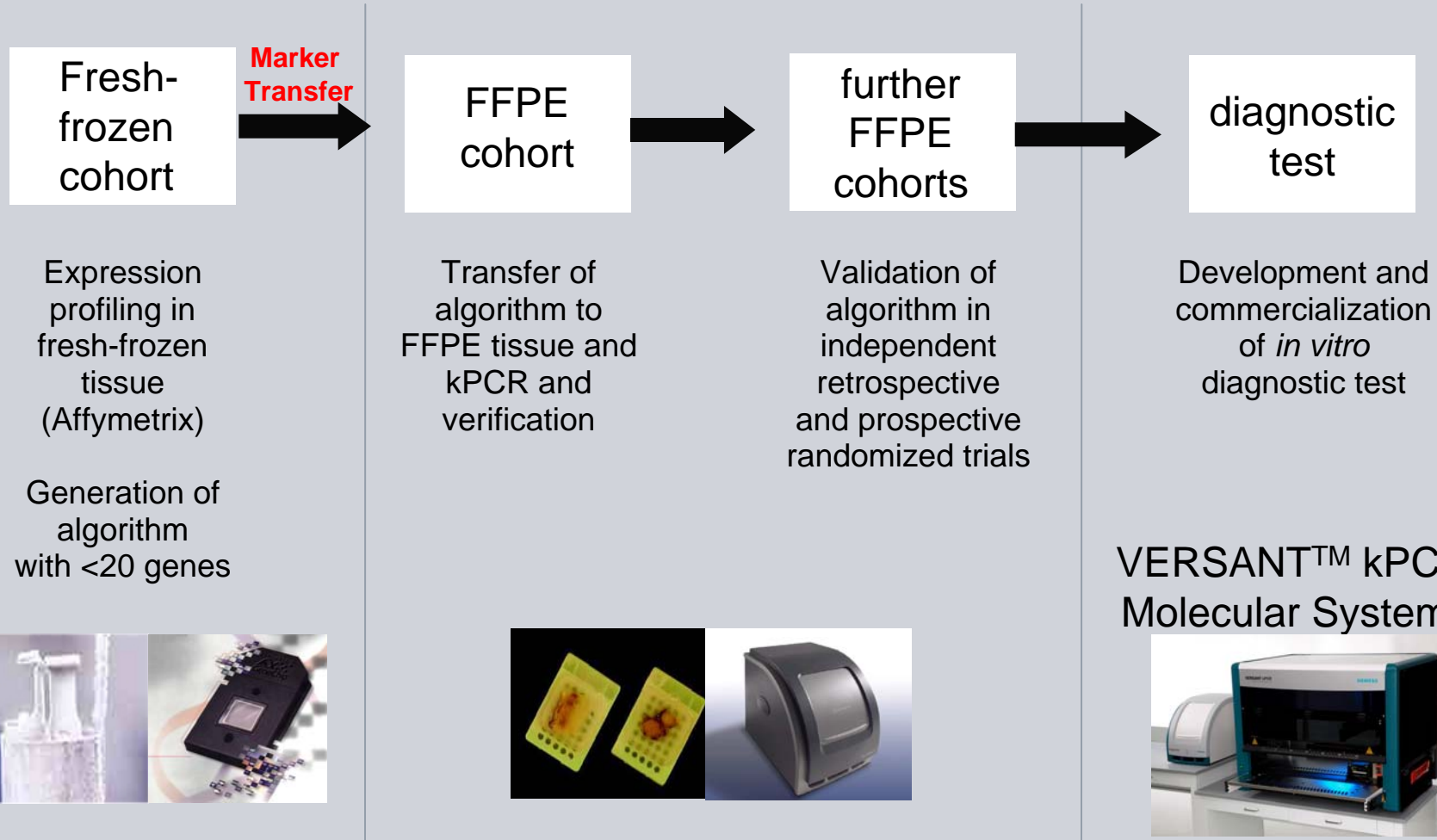

## Issues with Total RNA from FFPE Tissue

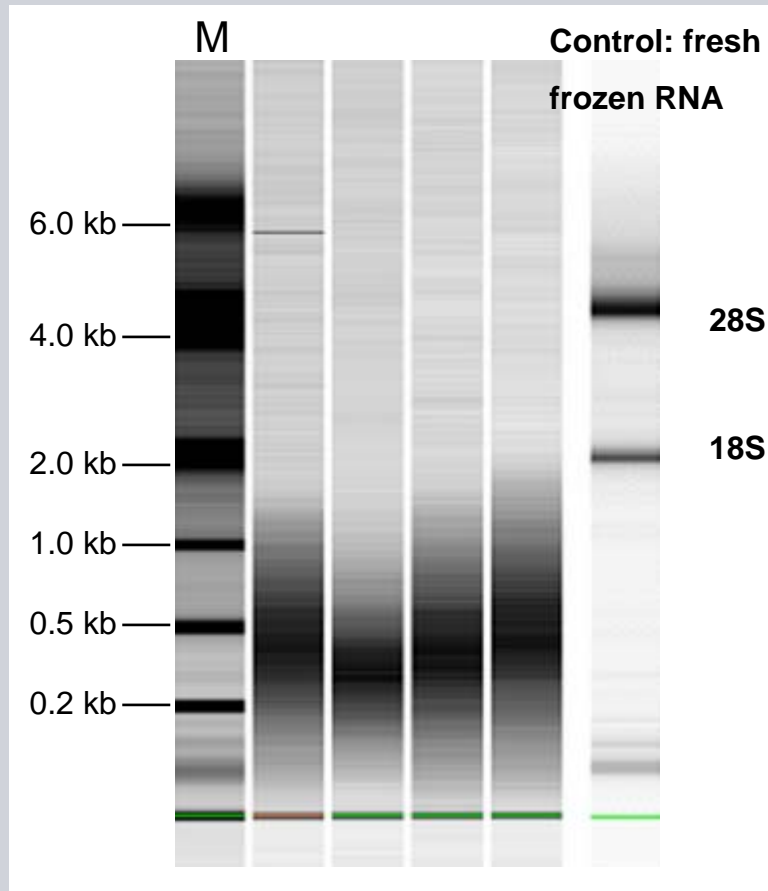

- RNA from fixed tissue on Bioanalyzer
- RNA is chemically modified and crosslinked to other nucleic acids and protein
- RNA is fragmented

← Mean size  
~ 200-500 bp

Fragmentation over storage time


## Open Questions for Accurate Expression Profiling in FFPE Tissue

**SIEMENS**


- No Standardized Isolation Technology for Nucleic Acids
- Different Fixation Conditions (Time in Fix / Time to Fix)
- Variability of Tumor Cell Content
- Block to Block Variability (Intra-tumor Heterogeneity)
- Correlation to Immunohistochemistry

## Siemens' „New“ Magnetic Particles

- „One particle chemistry“ for all applications in the field of nucleic acid isolation
- Coating of iron oxides with nanolayer of silica
- No change in morphology (spherical, evenly sized, superior active surface / volume ratio)

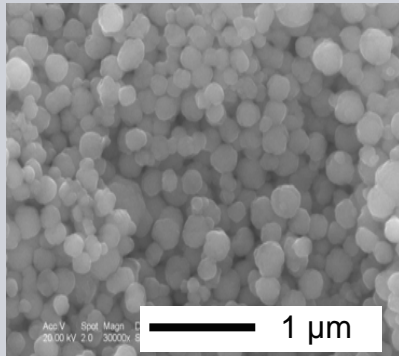

Siemens

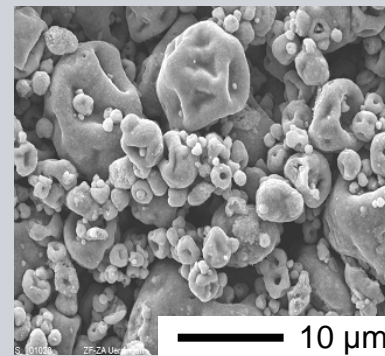


Others



# Siemens' Isolation Workflow for RNA (and DNA) from FFPE Tissue

SIEMENS

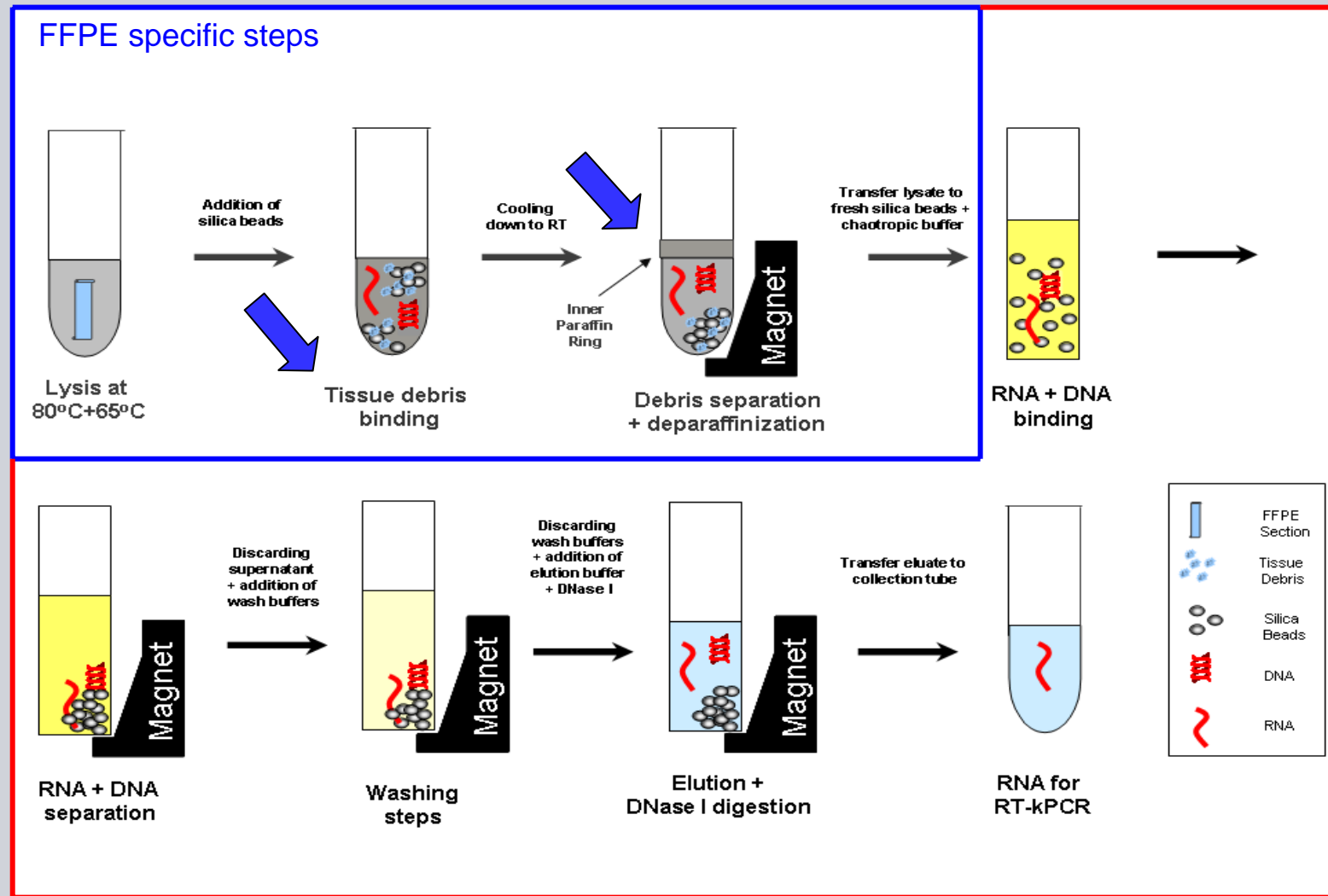

## Automation of DNA/RNA FFPE Isolation




### Adapted to VERSANT kPCR™ Platform

- Fully automated and high throughput (48 sections in 4 hours)
- 10 µm sections are loaded in plastic tubes on robot
- No manual deparaffinization with xylene/ethanol and centrifugation
- Standardization of lysis conditions by negative selection step for “tissue debris”
- Currently in development

**Performance (Yield of Total RNA from FFPE Sections)**

**Better Yield**

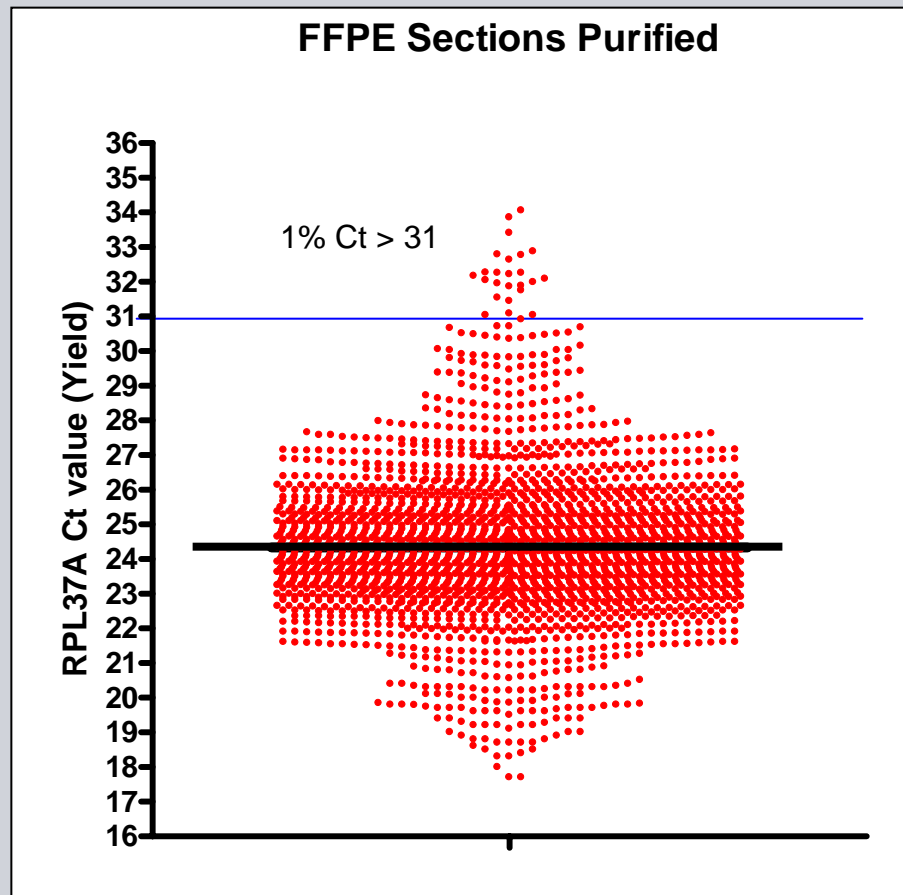Breast Cancer Tumor Blocks:  
3 month to 21 years of age

**n ≥ 2314**



Median RPL37A Ct = 24.5  
~ 1 µg / 10 µm section

**Success Rate: 100%**

**High Yield Rate: 99%**

Sample requirement for RT-kPCR: 2 ng/well ⇒ ~ 500 PCRs per slide

## Between Instruments Reproducibility of mRNA Expression

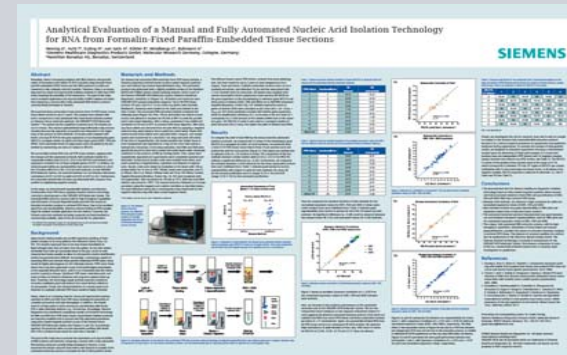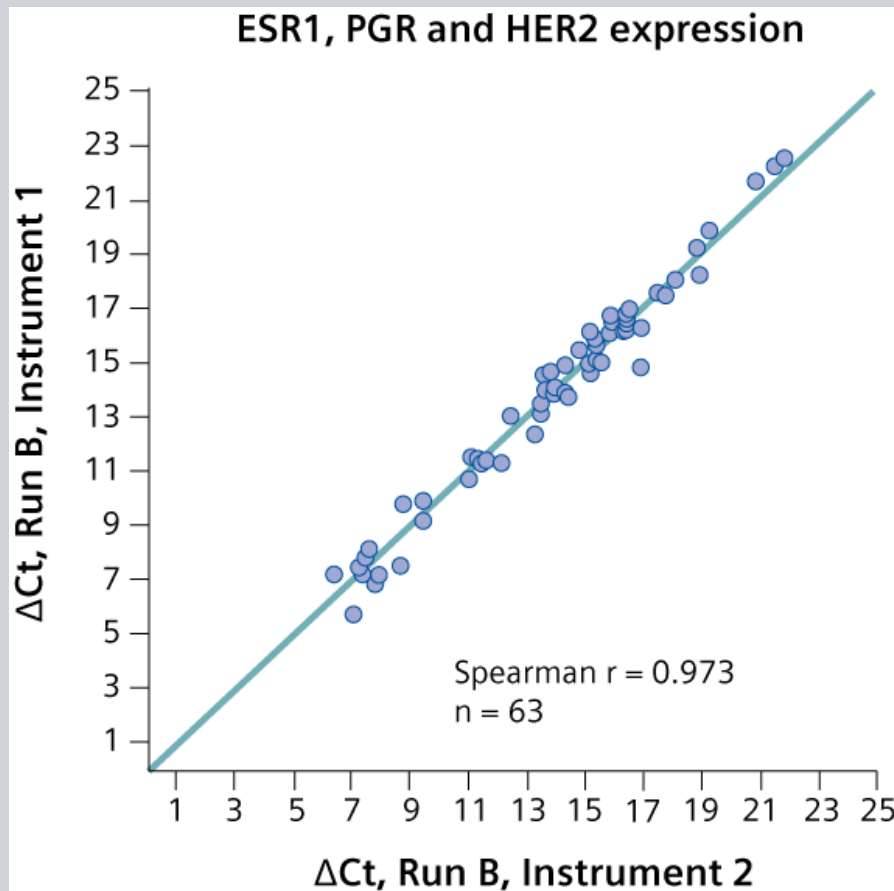


Poster AMP 2008

Robust and reproducible method without Carry Over of target RNA from high positive to neighbouring negative sections

**Benchmark to State-of-the-art Qiagen Kit's**

**Total and hands-on time per 90 samples**



Bohmann et al, 2009 (submitted)

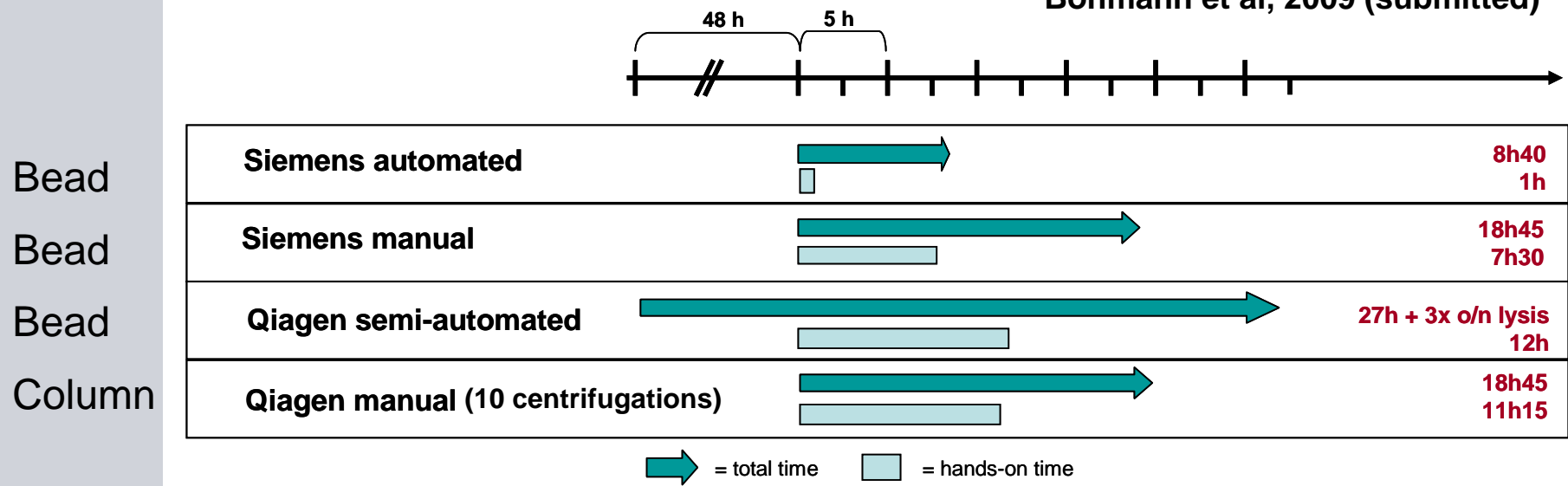

- Automated protocol superior with regards to total time and hands-on-time
- Automated protocol showed best section to section reproducibility



## Influence of Storage Time on Yield

Müller et al, 2009, submitted (cooperation with Institute of Pathology, Charite Berlin)  
 Extraction of RNA from 167 breast carcinomas stored 2 months up to 21 years

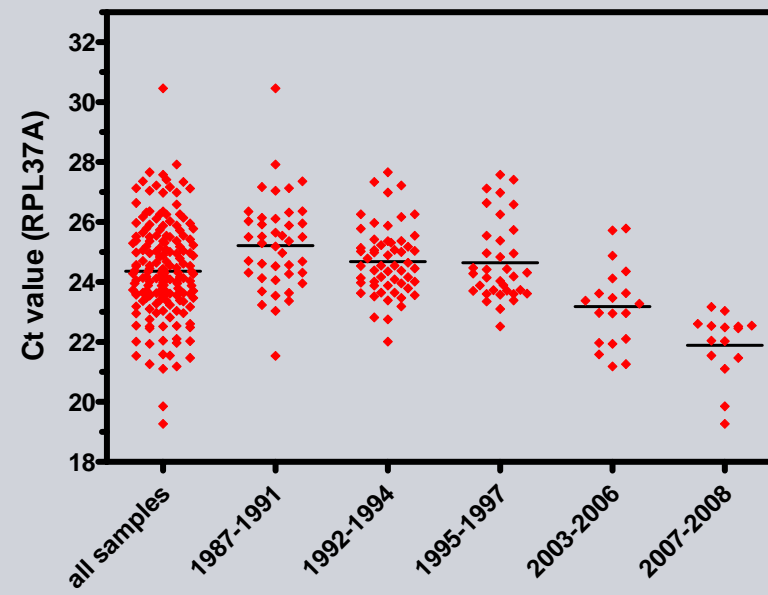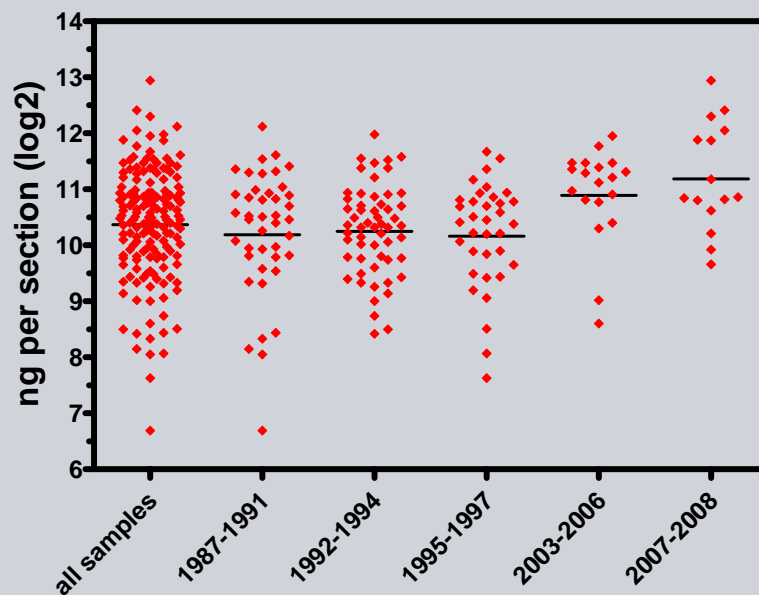



Yield of total RNA by Ribogreen or RPL37A Ct value significantly decreases with storage time

## Section to Section Reproducibility of Normalized Expression

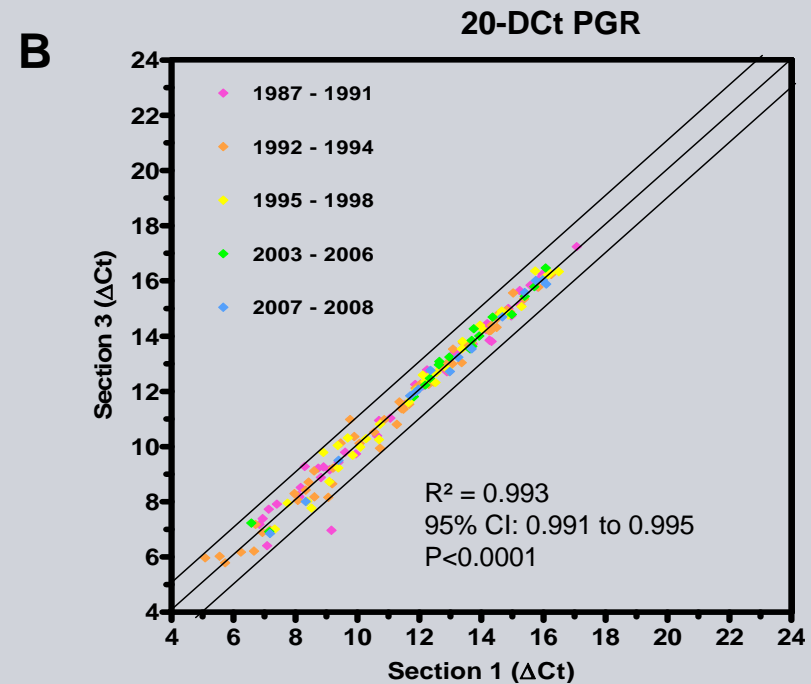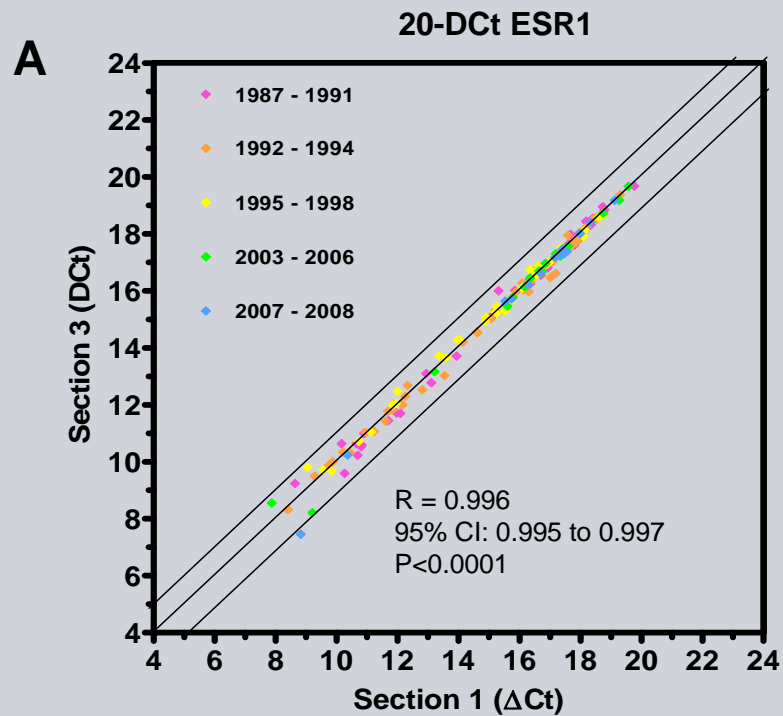

High reproducibility of quantitative RNA expression values ( $\Delta$ Ct for ESR, PGR and ERBB2) between consecutive sections

# High Concordance between Protein and RNA level

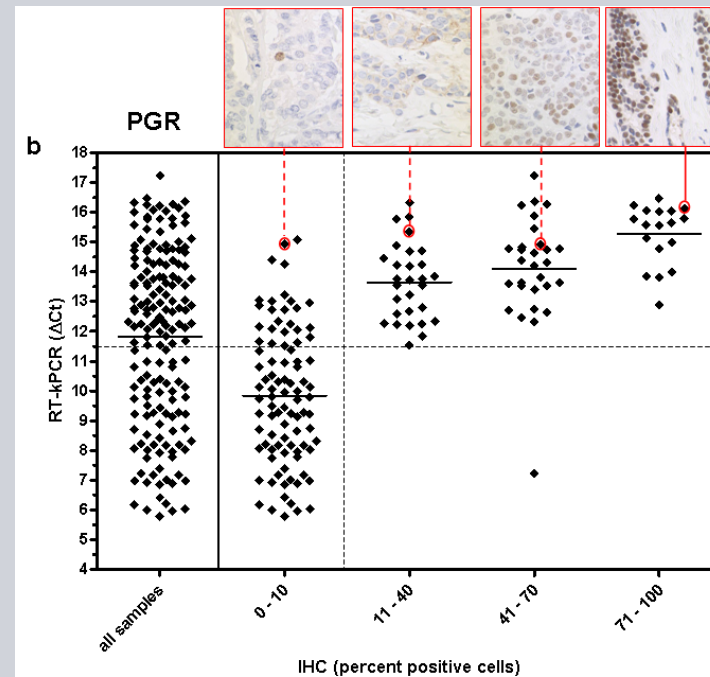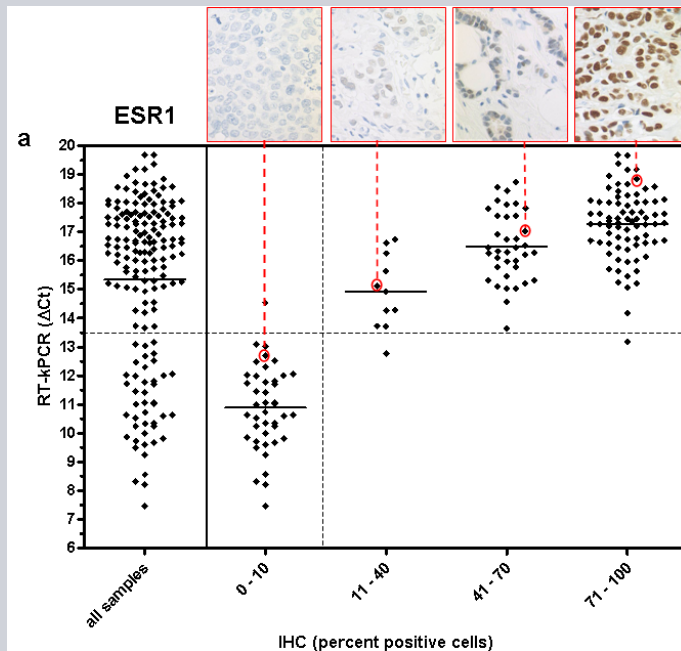

RT-PCR  
Cut-off

Overall Agreement	ESR1	98,2%	(Pos 98,4% / Neg 97,6%)
	PGR	84,4%	(Pos 98,7% / Neg 72,5%)

High concordance between quantitative RNA expression and semiquantitative immunohistochemistry data for ESR1 and PGR



## Different Time **in** Fixation Study

Cooperation with Prof. Porembe,  
Center for Histology, Cytology and Molecular Diagnostics, Trier

### 1 Breast Cancer Tumor

Portion 1: TIF 1h in Neutrally Bufferd Formalin (NBF)

Portion 2: TIF 20h in NBF (“routine”)

Portion 3: TIF 2 days in NBF

Portion 4: TIF 5 days in NBF


Portion 5: TIF 10 days in NBF

Portion 6: Fresh frozen within a few minutes (control)

Duplicates from each portion (independent isolations + detections)

10 genes detected with RT-kPCR and normalized to RPL37A

# Time in Fixation Normalized Expression Data

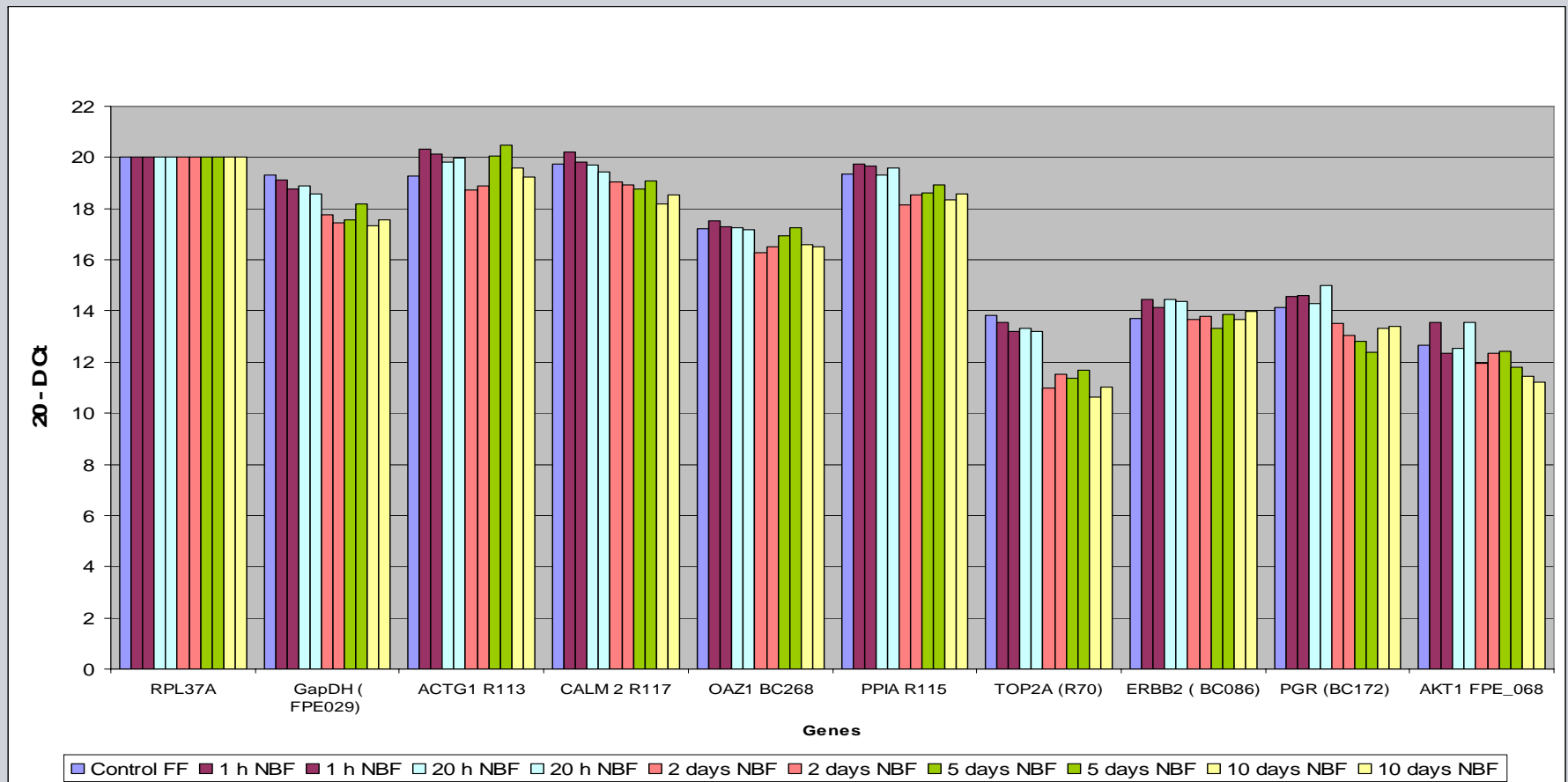

Even extreme fixation times allow accurate expression profiling in RNA from FFPE tissue

## Different Time **to** Fixation Study

1 Breast Cancer Tumor

Portion 1: TTF 10 min at 20°C (“routine”)

Portion 2: TTF 10 min at 37°C

Portion 3: TTF 1h at 4°C

Portion 4: TTF 1h at 37°C

Portion 5: TTF 12h at 4°C

Portion 6: TTF 12h at 20°C

Portion 7: TTF 12h at 37°C

Portion 8: TTF 24h at 20°C

Portion 9: TTF 24h at 37°C


Portion 10: FF within a few minutes (control)

**Portion 1-9 were fixed with  
neutrally buffered formalin  
(NBF) for 16-20h**

Duplicates from each portion (independent isolations + detections)

10 genes detected with RT-kPCR and normalized to RPL37A

## Different Time to Fixation Normalized Expression Data

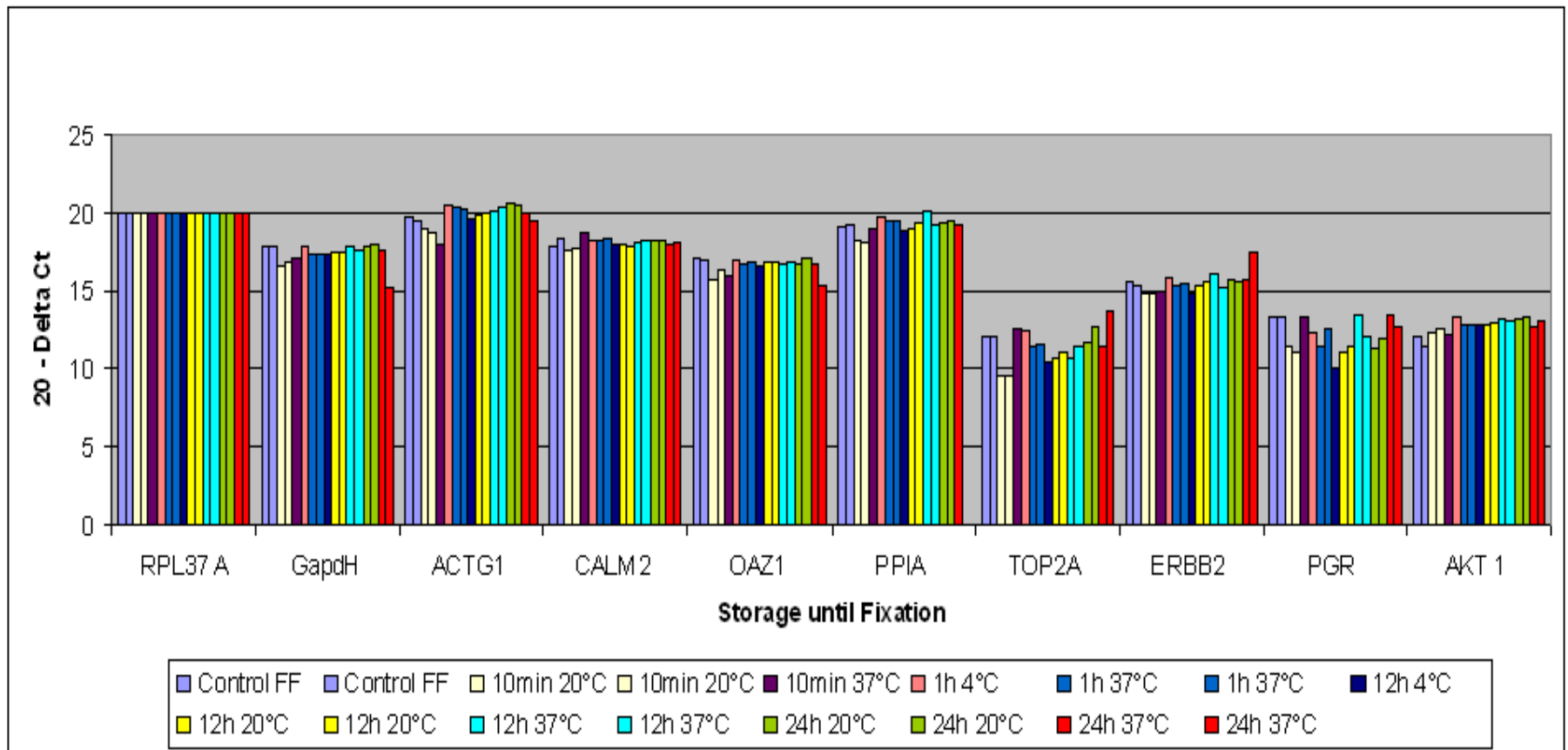


Even extreme time intervals from surgery to the fixative allow accurate expression profiling in RNA from FFPE tissue

## Tumor Content + Heterogeneity Study



Cooperation with T. Tramm + J. Overgaard,  
Department of Experimental Clinical Oncology, Aarhus University Hospital, Denmark

- Paraffinblocks from 22 pts. diagnosed with breast cancer 1991-1993
- Two independent FFPE blocks from one tumor
- 1 HE-section from each block
- H: whole section = 30 - 100 % tumor
- T+: trimmed section ( invasive component) = 90 - 100 % tumor
- Automated isolation with Siemens´ FFPE RNA method
- Genes detected with RT-kPCR: RPL37A, ESR, PGR, ERBB2

ASCO 2007

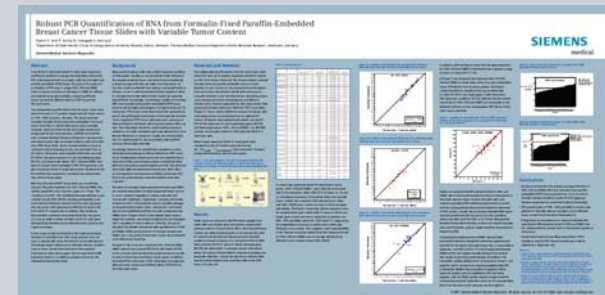


## Different Tumor Content Study

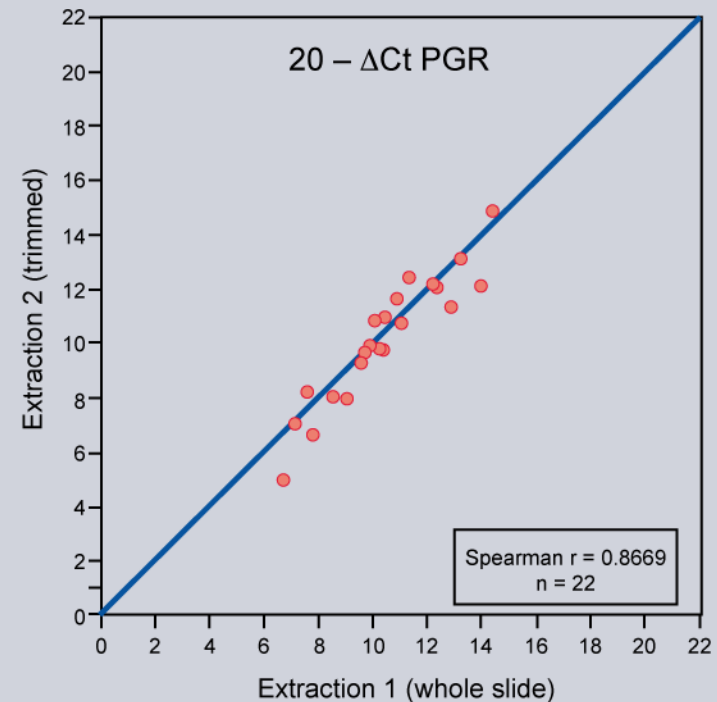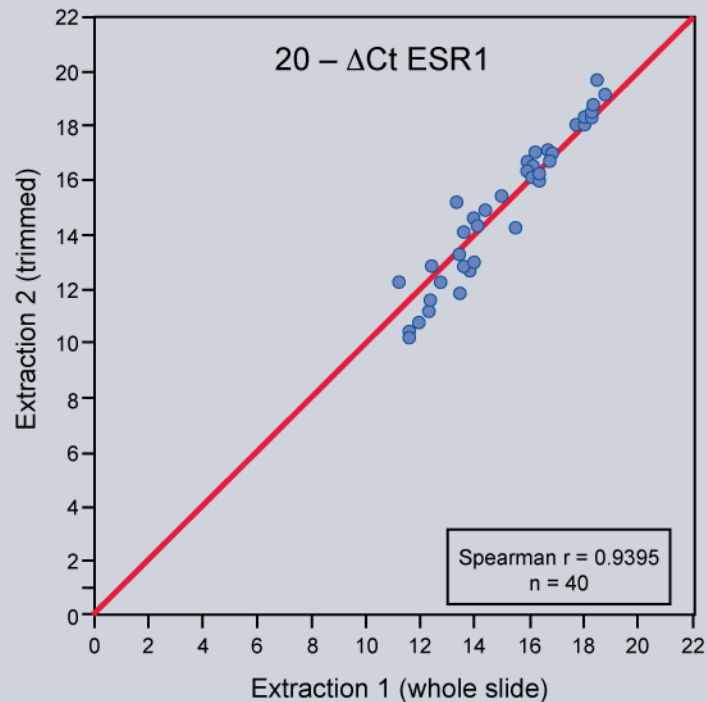

High reproducibility between independent isolations and different tumor content for ESR, PGR and ERBB2 RNA expression

## Intra-Tumor Heterogeneity

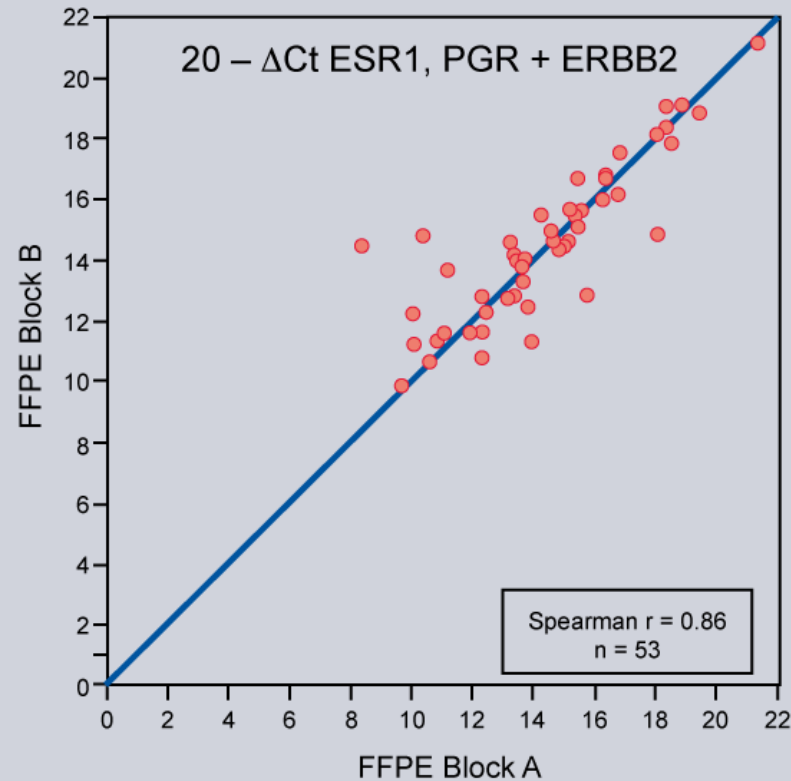

Good correlation between independent FFPE blocks from one patient for ESR, PGR and ERBB2 RNA expression

## Summary

- Siemens developed a unique nucleic acid isolation technology for RNA (and DNA) from FFPE tissue based on silica magnetic beads
  - High yield + Carry Over free
  - Fully automated (No deparaffinization / High throughput)
  - Reproducible and robust (under different clinical conditions)
- Technology is key for detection and quantification of molecular markers by expression profiling (PCR or arrays), sequencing and genotyping
- Contribution for implementation of robust and high-throughput testing of biomarkers in clinical routine and retrospective + prospective research studies



## Acknowledgements



K. Bohmann, R. Kronenwett, U. Stropp, C. Petry, K. Weber  
T. Acht, H. Euting, R. Köhler, M. Averdick, C. Windbergs,  
C. Roth, G. Altmann



University Hospital Aarhus

T. Tramm and J. Overgaard  
Department of Experimental Clinical Oncology; Aarhus



C. Denkert and B. Müller  
Institute of Pathology, Charite Berlin



K-L Schäfer, Uwe Rogel  
Institute of Pathology, Düsseldorf

Center for Histology,  
Cytology and Molecular  
Diagnostics

Ch. Poremba, Trier